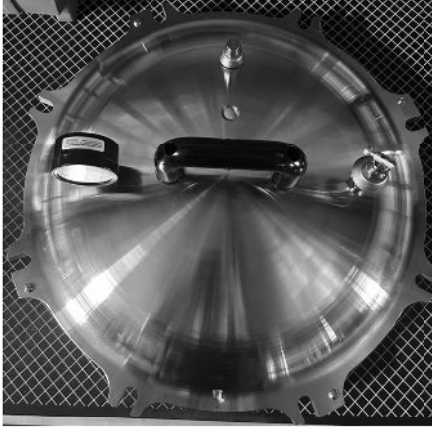
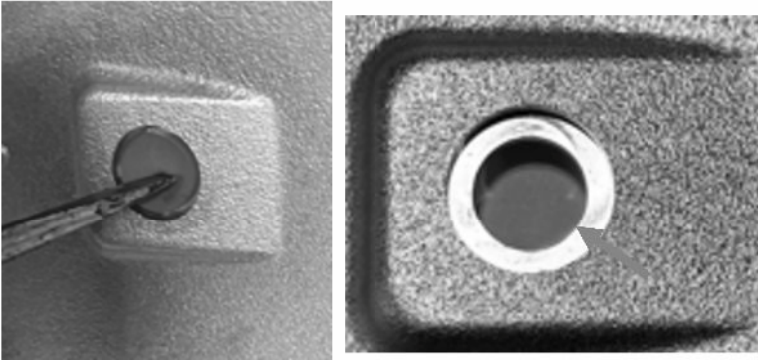


Component Replacement Instructions

Consumer Products
1931 South 14th Street
Manitowoc, WI 54220

Part Name	Overpressure Plug	Author	AD
Part Number	1010	Revision Date	02/01/2024
Instruction Details		Photos/ Images	
		Tools Needed: 1. Screwdriver	
Removal Instructions 1. Use screw driver to remove by pressing through from the back. WARNING  Careful not to damage the edges of the hole risk of not being able to maintain proper operating pressures.			
Installation Instructions 2. Install by starting on an angle from the back side then continue pressing through using your finger.			
3. Turn cover over and confirm that it is pressed through and that there are no edges of the plug folded under. WARNING  If not assembled correctly risk of the plug blowing out before operating pressure is achieved.			

Component
Replacement Instr

Consumer Products
1931 South 14th Street
Manitowoc, WI 54220



Part Name	Overpressure Plug	<i>AD</i>
Part Number	1010	02/01/2024
Instruction Details		